
PowerLab SP Owner's Guide

for PowerLab/4sp, PowerLab/8sp, and PowerLab/16sp models

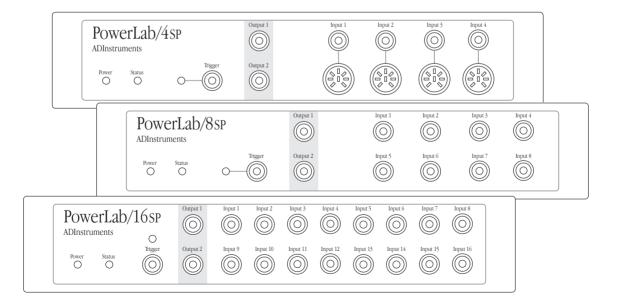

This document was, as far as possible, accurate at the time of printing. Changes may have been made to the software and hardware it describes since then, though: ADInstruments reserves the right to alter specifications as required. Late-breaking information may be supplied separately.

Trademarks of ADInstruments

MacLab, PowerChrom, and PowerLab are registered trademarks of ADInstruments Pty Ltd. E-Corder and the names of specific recording units, such as PowerLab/16SP, are trademarks of ADInstruments Pty Ltd. Chart, EChem, Histogram, Keeper, Peaks, Scope, Translate, UpdateMaker, and UpdateUser (application programs) are trademarks of ADInstruments Pty Ltd.

Other Trademarks

Apple, the Apple logo, Mac, Macintosh, and PowerBook are registered trademarks of Apple Computer, Inc. Power Macintosh, iMac, and System 7 are trademarks of Apple Computer, Inc.

PowerPC is a trademark of International Business Machines Corporation.

Pentium is a registered trademark of the Intel Corporation.

Windows and Windows 95 are registered trademarks of Microsoft Corporation. Windows NT is a trademark of Microsoft Corporation.

All other trademarks are the properties of their respective owners.

Product: PowerLab/4sp (ML750); PowerLab/8sp (ML785); PowerLab/16sp (ML795)


Hardware: Michael Macknight, Boris Schlensky, Tim Spencer, Graham Milliken, Indulus Kradzins.

Documentation by staff of ADInstruments Document Number: U-ML003S-OG-002E Copyright © September 2001 ADInstruments Pty Ltd Unit 6, 4 Gladstone Rd Castle Hill, NSW 2154 Australia

Web: http://www.adinstruments.com E-mail: support@adinstruments.com

All rights reserved. No part of this document may be reproduced by any means without the prior written permission of ADInstruments Pty Ltd.



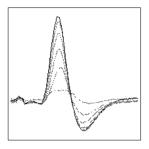

Contents

Contents iii

Safety Notes $\,\mathrm{v}$

1 Overview 1

How to Use this Guide 2 First, Check Your PowerLab! 2 The PowerLab System 3 Computer Requirements 3 The Application Programs 4 ADInstruments Front-Ends 5 The PowerLab 5 The Front Panel 5 The Back Panel 9

2 Setting Up 15

Self-Test 16 The USB Connection 17 Connecting the PowerLab Using USB 18 USB Connection Rules 19 The SCSI Connection 20 SCSI Connection Rules 20 Setting the SCSI ID Number 22 SCSI Termination Rules 23 Connecting the PowerLab: Macintosh 23 Connecting the PowerLab: Windows 24 SCSI Strikes Back 26

A Technical Aspects 27

How Does it Work? 27 The Analog Inputs 31 PowerLab Accuracy 33 The External Trigger 33 The Analog Output 34 Connections 35 USB Port 35 SCSI Port 35 I²C Expansion Port 36 Digital Input and Output Ports 37 Pod Connectors 37

B Troubleshooting 39

Problems: Macintosh 39 Problems: Windows 42

C Specifications 47

PowerLab/4sp Specifications 47 PowerLab/8sp Specifications 50 PowerLab/16sp Specifications 53

Glossary 57

Index 61

Licensing & Warranty Agreement 63



Product Intention

PowerLab systems have been designed only for use in teaching and research applications. They are not intended for clinical or critical lifecare use and should never be used for these purposes, nor for the prevention, diagnosis, curing, treatment, or alleviation of disease, injury, or handicap.

Applicable Safety Standards

When used with insulated transducers or ADInstruments isolated front-ends, PowerLab systems are safe for human connection. The ML132 Bio Amp, ML135 Dual Bio Amp, ML408 Dual Bio Amp/Stimulator, ML116 GSR Amp, ML117 BP Amp and ML180 frontends conform to international safety requirements. Specifically, these requirements are IEC601-1 and its addenda (Table S–1), and various harmonised standards worldwide (UL2601.1 in the USA, and AS/NZS3200.1 in Australia and New Zealand). In accordance with European standards they also comply with the electromagnetic compatibility requirements under EN60601-1, which encompasses the EMC directive. Further information is available on request.

Table S-1Applicable safety standardsmet by ADInstrumentsisolated front-ends

EC Standard No.	International Standard — Medical electrical equipment
IEC601-1:1988	General requirements for safety
IEC601-1-1:1992	Safety requirements for medical electrical systems
IEC601-1-2:1993	Electromagnetic compatibility

General Safety Instructions

- Since the PowerLab is used in conjunction with a computer (which has much higher leakage currents) the subject must be at least 1.83 metres (6 feet) away from any of the equipment, to avoid contact with the computer. The Bio Amp or BP Amp cable provides enough distance. Alternatively, an isolation transformer can be used to supply power to both the PowerLab and the computer.
- The PowerLab /4SP, PowerLab /8SP and PowerLab /16SP should always be fitted with the recommended fuses for safe operation. Specific information on fuse types and replacement is given in a later chapter.
- The PowerLab /4SP, PowerLab /8SP and PowerLab /16SP are classified as Class I medical equipment, which means that protection against electric shock in the event of a fault relies on a direct connection through the power cable to your building's earth conductor. The power cable supplied with your PowerLab provides the required ground connection to the power outlet. If your building does not have power outlet sockets with a good ground connection, then you may use the ground connection on the rear of the PowerLab to provide the equipotential connection to the building's earth conductor. A ground connection is an essential part of this equipment's safety. Never use the PowerLab without a ground connection.

Safety Symbol Explanation

Every ADInstruments device designed for connection to humans, including the Bio Amp, BP Amp, GSR Amp and Stimulus Isolator, carries one or more of three safety symbols, as shown in Figure S–1. These symbols appear next to those input and output connectors that can be directly connected to human subjects.

Figure S–1

Equipment safety symbols on ADInstruments devices designed for safe connection to humans



BF symbol: Bodyprotected equipment



documentation



CF symbol: Cardiacprotected equipment

The three symbols are:

- **BF (body protected) symbol**. This means that the input connectors are suitable for connection to humans provided there is no direct electrical connection to the heart.
- **Warning symbol**. The exclamation mark inside a triangle means that you should consult the supplied documentation (you're reading it) for operating information or cautionary and safety information before using the device.
- **CF (cardiac protected) symbol**. (Appears on some models of Bio Amp, and every BP Amp.) This means that the input connectors are suitable for connection to humans even when there is direct electrical connection to the heart.

Cleaning and Sterilisation

This system is not supplied with materials or components likely to need sterilisation. If you need to clean peripherals likely to be in contact with different subjects, they can be cold-sterilised with an appropriate sterilising agent. No part of the system can be autoclaved.

Storage

It is recommended that the hardware be kept below 40 °C and above 0 °C in a moisture-free or low-humidity environment when not in use (storage only). Electronic components are susceptible to corrosive substances and atmospheres, so it is also advisable to keep the system away from laboratory chemicals.

Preventative Inspection and Maintenance

Both PowerLab systems and ADInstruments front-ends are maintenance free and do not require periodic calibration or adjustment to ensure medical safety. Internal diagnostic software performs system checks during power up and will report errors if a significant problem is found.


There is no need to open the instrument for inspection or maintenance, and doing so within the warranty period will void the warranty.

If you so wish, your PowerLab system can be periodically checked for basic medical safety by using an appropriate medical safety testing device. Tests such as earth leakage, earth bond, insulation resistance, patient leakage and auxiliary currents and power cable integrity can all be performed on the PowerLab system without having to remove the covers. You should follow the instructions for the testing device if performing such tests.

If the PowerLab system is found not to comply with such testing you should contact your PowerLab dealer to arrange for the equipment to be checked and serviced. You should not attempt to service the device yourself.

1

CHAPTER ONE Overview

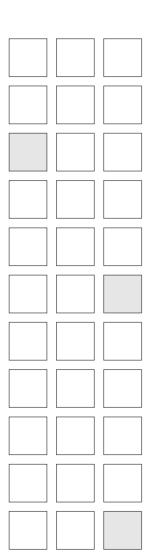

Your PowerLab recording unit, together with a range of specialised application programs, provides a versatile data recording and analysis system when used with a Windows or Macintosh computer. This chapter provides an overview of the PowerLab system and describes the basic features, connectors, and indicators of the SP series PowerLabs: the PowerLab/4SP, PowerLab/8SP, and PowerLab/16SP.

How to Use this Guide

This owner's guide describes how to set up and begin using your PowerLab recording unit. The chapters provide an overview of the PowerLab system (the combined software and hardware package), and a more detailed look at the features of your recording unit and its connection to your computer. The appendixes provide technical information about the recording unit, and solutions to problems. At the end of this guide is a glossary of hardware terms, an index, and warranty information.

The specifications and diagrams included in the appendixes are there to help the more technically minded to understand what the PowerLab can and cannot do, but this is not a service manual: only an authorised ADInstruments distributor should attempt repairs. If you modify the recording unit yourself, you void any rights you have under warranty.


The user's guides for the Chart and Scope application programs provide detailed information on the software side of the PowerLab system and its uses in acquiring, storing, and analysing data. Read them after you have connected the PowerLab to your computer.

First, Check Your PowerLab!

Please do not attempt to connect the PowerLab to a power outlet or computer or turn it on until you have read the first two chapters of this owner's guide, and have checked it as described below.

- 1. Check that all items in the accompanying packing list are included in the box.
- 2. Check that there are no obvious signs of external damage to the PowerLab.
- 3. Check that there are no obvious signs of internal damage, such as rattling. Pick the PowerLab up, tilt it gently from side to side, and listen for anything that appears to be loose.
- 4. Check that the correct voltage for your country is shown on the back of the unit. Your PowerLab should be delivered with the

Figure 1–1 The voltage rating on the back panel of the PowerLab

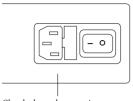

Check the voltage rating beneath the power switch

appropriate mains power voltage set, at either 100–120 volts or 220–240 volts. The setting is indicated at the right of the back panel of the unit, beneath the power switch.

If anything is missing, or the PowerLab seems to be damaged in any way, or if no voltage, or the wrong voltage, is shown on the back panel of the unit, contact your authorised ADInstruments distributor immediately, and describe the problem. Arrangements can be made to replace or repair the PowerLab. (Up-to-date contact addresses are available in the software — see Appendix B in one of the software user's guides for details.)

The PowerLab System

The PowerLab system is an integrated system of hardware and software designed to record, display, and analyse experimental data. The hardware consists of the PowerLab recording unit and possible ancillary devices (front-ends); the software consists of the application programs (such as Chart and Scope) that run on the computer to which the PowerLab is connected.

Your SP series PowerLab has considerable computing power of its own and performs many tasks that are necessary during data recording. Once the PowerLab transfers the data to the computer, the data are available for display, manipulation, printing, storage, and retrieval. The PowerLab/4SP has four inputs for recording external signals, the PowerLab/8SP has eight, and the PowerLab/16SP has 16.

Computer Requirements

Computer requirements for both Windows and Macintosh depend on whether you connect your /4SP, /8SP or /16SP PowerLab unit to your computer using a SCSI or a USB interface. SCSI allows faster transfer rates, but is less straightforward to use.

Windows

- A computer with a Pentium processor or better
- 32 MB free RAM (48 MB for Windows NT or 2000)

- Windows 95, 98, Me, NT 4.0, 2000 or later Windows operating system
- Additional requirements for USB: A USB interface on your computer (this may be added via a card if necessary); Microsoft Windows 98, Me, 2000, or later (Windows 95 and NT will not work with USB).
- Additional requirements for SCSI: An appropriate SCSI card and cable (we recommend that you use a card and cable supplied by ADInstruments — others may be suitable, but you should seek technical advice if in doubt); a software component called ASPI (ASPI will already be on your computer unless you are using Windows NT or 2000, in which case you can install it using the installer located on the PowerLab Installer CD).

Macintosh

These requirements assume you are using the latest version of Chart. An older version of Chart is available for use with older Macintosh computers.

- A Power Macintosh computer (or computer with a PowerPC microprocessor)
- 32 MB free RAM (64 MB recommended)
- Mac OS 8.6 or later (Mac OS 9.0 or later recommended)
- Additional requirement for USB: A USB interface on your computer (this may be added via a card if necessary)
- Additional requirement for SCSI: Either built-in SCSI or an appropriate SCSI card and cable (we recommend that you use a card and cable supplied by ADInstruments others may be suitable, but you should seek technical advice if in doubt).

The Application Programs

Two main application programs and their documentation are provided with each PowerLab. Scope emulates a two-channel storage oscilloscope. Chart emulates a multi-channel chart recorder (up to 16 channels, depending on the PowerLab model). Both provide many other powerful features in addition, including computed functions, triggering options, software-controlled stimulus generation, and automated recording and analysis. Their user's guides describe them in full; the Getting Started with PowerLab manual gives summaries. You must install the software to use your PowerLab. Installation instructions are included in the user's guides for the software.

ADInstruments Front-Ends

Front-ends are optional ancillary devices that can be connected to the PowerLab to extend the system's capabilities. They afford additional signal conditioning and other features, and extend the types of experiments you can conduct and the data you can record. For instance, the Bio Amp is available if you need to perform electrically isolated measurements of biological signals (PowerLabs should *never* be connected to humans directly, unless it is through connections marked as safe). A PowerLab can have as many front-ends connected to it as it has analog inputs, roughly speaking. Front-ends are easily added to or transferred between PowerLabs. They are automatically recognised by the PowerLab system and seamlessly integrated into its programs, operating under full software control.

Your ADInstruments distributor will have full information on available front-ends and their uses and specifications, and will be happy to discuss your requirements.


The PowerLab

It is a good idea to get familiar with some of the external features of your PowerLab before connecting it to a power source. The rest of this chapter discusses the different features, connectors, and indicators of the PowerLab/4SP, PowerLab/8SP, and PowerLab/16SP.


The Front Panel

The front panel of your PowerLab provides most of the connectors for interfacing with external signals, and indicators for various functions. This section describes each of the front panel features.


The PowerLab/4SP has three indicators at the left of the front panel, one BNC connector for the external trigger, two BNC connectors for output, and four BNC connectors (marked Input 1 to Input 4), with four alternative Pod (DIN) connectors, for recording external signals.

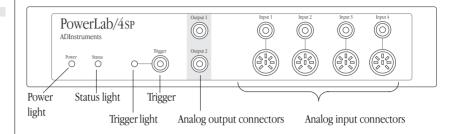

The PowerLab/8SP has three indicators at the left of the front panel, one BNC connector for the external trigger, two BNC connectors for output, and eight BNC connectors (marked Input 1 to Input 8), for recording external signals.

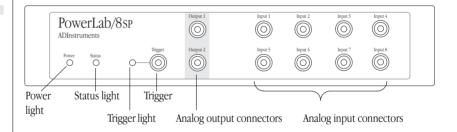


The PowerLab/16SP has three indicators at the left of the front panel, one BNC connector for the external trigger, two BNC connectors for output, and 16 BNC connectors (marked Input 1 to Input 16), for recording external signals.

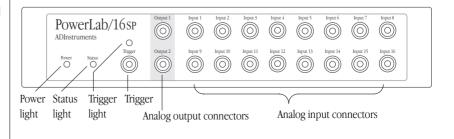





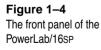

Power and Status Indicators

The Power indicator on the front panel is a green light, which simply shows that the PowerLab is getting power. The indicator next to it, the Status indicator, provides a visual indication of what the PowerLab is doing, and will flash different patterns and colours depending on the state of the PowerLab.

Status Indicator	Meaning
Off	Idle and not yet initialised by the software.
Yellow flashes	When the PowerLab is powered up, the indicator flashes yellow a number of times equal to the SCSI ID. For example, 4 yellow flashes would indicate a SCSI ID number of 4.
Green constant	Idle, initialised, and waiting for a command from the computer.
Yellow constant	Sampling, or communicating with the computer.
Yellow flash then one or more red flashes	This indicates that the PowerLab has detected an internal fault during the power-up test. The number of red flashes after the yellow flash indicates the type of error that has been detected. This display will keep repeating until the power to the PowerLab is turned off.

Analog Inputs

The analog inputs used to record external signals can handle signals from ± 10 V down to the microvolt (μ V) range without the need for additional external amplification. Each PowerLab analog input has an independently programmable gain amplifier with its own filtering, and AC/DC coupling. You can set up each input independently to suit your requirements using the software. Note that applying more than ± 15 V to the analog inputs can damage the circuitry.

The PowerLab/4sP has four independent analog inputs (marked Input 1 to Input 4), each with alternative connectors. The top BNC connector can be used for single-ended input, and the 8-pin DIN Pod connector below can be used for either single-ended or differential input. Single-ended inputs record the difference between signal and ground, and differential ones record the difference between positive and negative input signals. Do not attempt to record from both the BNC and Pod connectors for an input at the same time, or the signals will compete.

 Table 1–1

 Status indicator states (don't

worry too much about these)

▲ Caution

PowerLab inputs and outputs are not electrically isolated. Human subjects must not be connected directly to the PowerLab. If such measurements are to be made, an isolated front-end must be used, such as the Bio Amp. Pod connectors allow the connection of ADInstruments Pods small, low-cost units that provide alternatives to front-ends for specific tasks, for use with precalibrated transducers and so on. (Pods are supported by Scope v3.6.3 and later, Chart for Macintosh v3.6.3 and later, and Chart for Windows v3.4.7 and later.) The Pod connectors on the PowerLab/4SP do not handle transducers directly unless the transducers are so labelled (unsuitable transducers will give a very weak signal). Transducers designed for direct connection can be provided with power and control, since the Pod connectors provide some functions of the I²C output as well as alternative analog inputs to the BNC connectors.

The PowerLab/8sP has eight independent analog inputs (marked Input 1 to Input 8). All inputs are single-ended (with the signal referenced to ground), and have only positive connectors.

The PowerLab/16SP has 16 independent analog inputs (marked Input 1 to Input 16). All inputs are single-ended (with the signal referenced to ground), and have only positive connectors.

Analog Output

The PowerLab can generate a stimulus voltage through its analog output sockets (marked Output 1 and Output 2), giving positive, negative, differential, or independent stimuli, depending on the sockets used and the software settings. By default, the outputs are used for complementary (differential) stimulation, where Output 1 is positive and Output 2 is negative. When Output 1 is used, a positive stimulus voltage (set up in the software) gives a positive voltage output, and a negative voltage a negative one. When Output 2 is used, the voltage outputs are inverted. When both output 2 is used, the stimulus is the difference between the voltages at the positive and negative outputs: you could generate up to a 20-volt pulse, given a ± 10 V stimulus.


The outputs can be used for independent stimulation in Scope and the Macintosh version of Chart, but not in the current version of Chart for Windows (v3.4). The details are discussed in the software user's guides.

Trigger

The external trigger connector allows you to use a digital signal level to synchronise recording to an external event. This input can handle voltages of up to ± 12 V. The threshold voltage (the voltage above which the trigger circuit activates) is 2.9 volts for the SP series PowerLabs. When the trigger threshold is crossed, the indicator beside the external trigger connector will glow yellow. The external trigger is described in more detail in Appendix A, and the software documentation covers its practical use in normal recording.

The Back Panel

The PowerLab back panel provides the sockets to connect the PowerLab to the computer, front-ends, the power outlet, and so on. This section describes each of the back panel features. The back panels are the same for all the SP series PowerLabs.

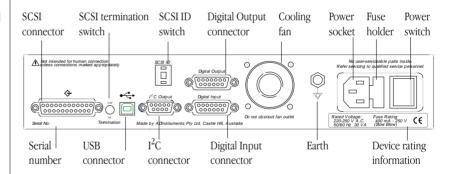

The USB Port

The PowerLab can connect to your computer using a USB (universal serial bus) port, which provides an alternative to the SCSI connection. This works with USB 1.1-compliant computers with USB connectors or a USB card. USB needs a Power Macintosh with Mac OS 8.5 or later, or a PC with Windows 98 or later.

A USB connection is a bit slower than SCSI for data transfer, and so may limit sampling rates (in Chart) or increase delay times between sweeps (in Scope), but it is less fussy. You can safely turn on or off, or disconnect or reconnect, a USB-connected PowerLab while the

Figure 1–5 The back panel of the PowerLab/4SP, PowerLab/8SP, and PowerLab/16SP computer remains on (this is *not* the case with SCSI). The application program (Chart or Scope) should not be on while this occurs, though. Please read the details on USB in Chapter 2 of this guide before connecting your PowerLab to your computer using USB.

The SCSI Port

The PowerLab can connect to your computer using a SCSI port, which provides an alternative to the USB connection. (SCSI is an acronym for small computer system interface, and is pronounced 'scuzzy' — which shows that engineers have no taste in acronyms). SCSI provides fast data transfer, but requires that specific rules be followed in order for it to work correctly: it is the most complex aspect of setting up a PowerLab system. Please read the details on SCSI in Chapter 2 of this guide before connecting your PowerLab to your computer using SCSI.

Each PowerLab has a single 25-pin connector for connecting SCSI cables, so it has to be used at the end of a SCSI chain. It is terminated internally, but the termination can be switched on and off using the SCSI Termination switch beside the SCSI connector (it should be on).

The SCSI ID Switch

The SCSI ID switch is there if you need to change the SCSI ID number of your PowerLab. The PowerLab is factory-set to an identification number of 4. The SCSI ID can be changed simply by pressing the control buttons while the PowerLab is turned off. The new ID takes effect when the PowerLab starts up again. Only some numbers are valid: they depend on the type of computer, whether it has built-in SCSI or a SCSI card, and the number of built-in and peripheral SCSI devices in the SCSI chain. This is discussed in the next chapter.

I²C Output

The I²C output is a special port designed to connect to front-ends made by ADInstruments. It supplies power and communications. A PowerLab can have as many front-ends connected to it as it has analog inputs, roughly speaking. *Note*. You should not attempt to run other external devices from the I²C port: it is designed for use only with ADInstruments front-ends. Only 50 mA maximum current can

be provided through this bus, so it should not be used for third-party devices drawing more current.

Digital Input and Output Ports

The digital input and output connectors let you monitor and control external devices respectively with the PowerLab. These recording automation and control features are available in the Macintosh version of Chart, but not in the current version of Chart for Windows (v3.4), or in Scope.

The digital input monitors state changes: you can have a predefined comment automatically inserted during recording when a digital input changes to a particular state. The eight lines of the connector allow monitoring of up to eight devices. The digital output can turn on and off external devices, for example, pumps, relays, and indicator lights, or can signal to some other device. The eight lines of the connector allow control of up to eight devices.

Technical details of the digital input and output connectors are given later in this guide: please refer to the descriptions before attempting to connect equipment to them. More information on the use of digital inputs and outputs is given in the user's guide for the software.


Ground Connection

A special earthing (grounding) stud has been provided on the rear of the PowerLab. This is an equipotential bonding connection post compatible with the DIN 42801 standard. If the ground connector post is used, the power cord ground connection should not be used: a power cord with no ground connection should be used instead.

The ground connector post on the rear panel is used as a primary earth connection (equipotential connection point) in situations that require this type of connection, or if there is no ground provided via the power cord. Safety standards in laboratories and similar environments may require additional grounding protection when connecting equipment to human subjects. In such cases, an equipotential connection may be used for all linked equipment to prevent ground loops, and power cords with no ground connections should be available.


Power Connections

The power switch on the back right of the PowerLab turns the PowerLab on and off; the 3-pin IEC power socket is used to connect your PowerLab to a 3-pin earthed (grounded) power cable.

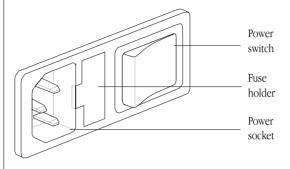

Fuse Replacement

The fuse holder is located beside the power socket, and can only be opened when the power cable is removed. To check the fuses, remove the power cable from the PowerLab, and lever the fuse holder out using the tab at its left, which is recessed in the power socket. The fuse holder contains two 'slow-blow' fuses, which protect your PowerLab from power surges.

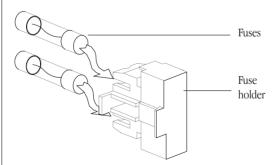

Replacement fuses must be of the correct type, as specified below for 100–120 volts or 220–240 volts operation:

100–120 V 1 Amp (5 × 20 mm) slow blow (voltage rating 120 or 250 V)

Figure 1–6 The power switch, fuse holder, and power socket

Figure 1–7

To change fuses in the PowerLab, remove blown ones from the fuse holder and slide new ones into place in the empty sockets • 220–240 V 400 mA (5 × 20 mm) slow blow (voltage rating 250 V only)

The correct voltage for your country is shown on the back of the PowerLab, by the power switch (see Figure 1–1, page 3).


Cooling Fan

The SP series PowerLabs can generate a fair amount of heat, so all of them are fitted with a cooling fan. The cooling fan is visible in the rear panel of the PowerLab. It maintains sufficient airflow in the enclosure to keep the circuitry within operating tolerances. The fan outlet must be kept unobstructed, and far enough from walls or other devices (at least three inches, 8 cm) so as to not impede ventilation.

It is also important to make sure that the air inlets underneath the PowerLab are unobstructed. Blocking the grilles may result in the PowerLab heating up enough internally to affect the performance of the system and reduce reliability.

2

снартек тwо Setting Up

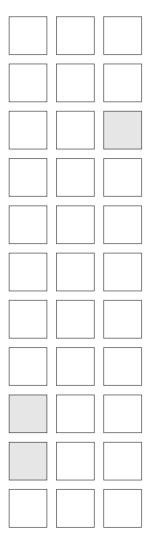

This chapter starts with the PowerLab's internal self-test, then looks at the USB connection. It also looks at the SCSI connection in some detail, including connection and termination rules, and setting the SCSI ID number. It discusses how to connect up your PowerLab to a Windows or Macintosh computer.

Self-Test

Now that you are familiar with some of the features of your PowerLab, you should check that it is working properly before you connect it to your computer.

The PowerLab performs a diagnostic self-test each time it is switched on, whether or not it is connected to a computer. To test that your PowerLab is functioning properly when you turn it on, follow the instructions below, and observe the effects.

- 1. Connect the PowerLab to a power outlet using the power cable that came with your unit. Turn the power on at the wall.
- 2. Turn on the power switch located on the rear of the unit.

The Power indicator on the front panel should glow green while the PowerLab is on (see Figure 1–2, Figure 1–3, and Figure 1–4).

The Status indicator should flash yellow a number of times equal to the SCSI ID of the PowerLab (see Table 1–1). If the internal diagnostic check finds no problems, the Status indicator turns off (showing that the PowerLab is idle; it only goes green when the software is running).

If the indicators perform as described above, then your PowerLab has successfully performed its internal self-test, and it can be safely connected to your computer. If your PowerLab would not successfully complete its self-test, something is wrong. If the Power indicator does not glow green when the power switch is turned on, then either the fuse has blown or there is a problem with the power cable or PowerLab itself.


If an error is detected during the self-test, the Status indicator will flash yellow and then a certain number of red flashes depending on the fault. Flashing lights are used when the problem is one that prevents the PowerLab communicating with the computer (such as a hardware fault), and so cannot be shown in the software. The error pattern will continue to be displayed, allowing you to take note of the type of error that occurred. The number of red flashes indicates the probable problem area. If the pattern is simply an alternating red flash and yellow flash, then the SCSI ID has been set to the numbers 0 or 7, which cannot be used for an SP series PowerLab. Switch off your PowerLab, change the SCSI ID for the PowerLab using the control on its back panel, wait about five seconds, and then switch it back on.


If the error pattern is not simple alternating red and yellow flashes, then something may be wrong with the hardware itself. The pattern of red flashes indicates the likely problem area to ADInstruments technicians. Take note of pattern, switch off your PowerLab, wait about five seconds, and then switch it back on. If it still continues to flash the pattern, then the PowerLab needs repair. Contact your authorised ADInstruments distributor as soon as possible. Do not attempt to repair the PowerLab yourself. If your PowerLab has successfully performed its internal self-test, read on to find out how to connect it to your computer.

The USB Connection

This section is essential reading before trying to use your PowerLab with USB. The PowerLab is fitted with a USB (universal serial bus) port, which provides an alternative to the SCSI connection. This works with USB 1.1-compliant computers with USB connectors or a USB card. On a PC, USB requires Windows 98, Me, 2000 or later (95 or NT 4 won't do). On a Macintosh, USB needs a Power Macintosh with Mac OS 8.5 or later (but note that the latest PowerLab software requires Mac OS 8.6).

USB devices are connected in a tree-like structure, with the host, the computer, forming the root of the tree, and hubs allowing multiple devices to connect to them, like the branches of the tree.

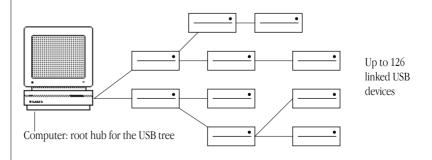

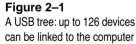


A hub is simply a device that lets you extend the USB tree. It connects to a USB device such as the computer, and multiple USB devices (including other hubs) can connect to it in turn. It provides power for those attached devices that need it. Up to five hubs can be connected between the computer and the furthest USB device. As many as 127 devices can be connected on one tree, counting the computer. The cable between any USB devices (including hubs) must allow a signal to be transmitted in a certain time; in practical terms this means it must be no more than 5 metres (16 feet) in length. Thus, no device can be more than 30 metres (98 feet) from the computer.

USB is more tolerant than SCSI. It lets one plug in and remove devices while the computer is on. It re-enumerates (provides addresses for) devices as they are connected and disconnected, rather than requiring fixed ID numbers. It has only one pair of connector types, unlike SCSI, and does not need termination as SCSI does.

USB still has some restrictions. A USB connection is a bit slower than SCSI for data transfer, and so may limit sampling rates (in Chart) or increase delay times between sweeps (in Scope). Since all the devices on a USB tree have a common communication pathway, if you have many devices with a lot of information to transfer, they compete for time. Using a video camera and a scanner at the same time as a PowerLab may limit the sampling rates considerably.


USB has two sorts of cables, which it should not be possible to interchange. Low-speed cables are cheaper, and suitable for mice and so on. They are either captive (one end is permanently fixed to the device) or use a custom connection, not a standard USB one. All detachable USB cables are high speed, with fully shielded twistedpair cables, and standard USB connections: a narrow rectangular A plug at one end and a squarer B plug with a bevelled top at the other. All devices that can act as hubs have high-speed connections. The PowerLab needs a high-speed connection, and is provided with a suitable cable.

Connecting the PowerLab Using USB

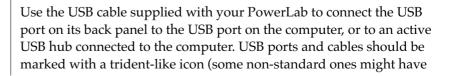



the letters 'USB' instead). USB cables are directional, and can only be connected one way: the narrow rectangular A plug connects to a hub (including the computer), and the squarer B plug with the bevelled top connects to a USB device, such as the PowerLab. Connecting Macintosh or Windows computers should be much the same in this case: just look for the icons. If the connection is to a USB card installed in the computer, there might not be any icons on the card. The narrow rectangular sockets for the USB A plug are distinctive, though. Any USB icons on the plugs should be on top.

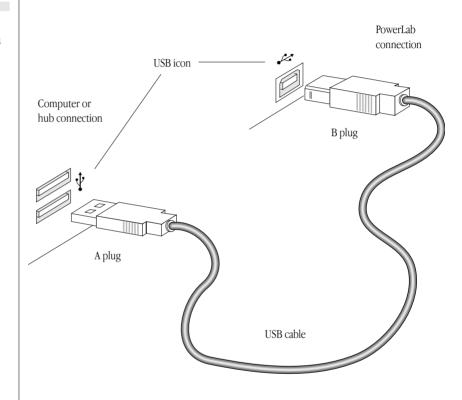

USB Connection Rules

 Cable length should be less than 5 metres (16 feet) between devices. Devices include USB hubs. There are a maximum of five hubs on any branch, so the maximum distance between the computer and the Powerlab is 30 metres (98 feet). Do not use extension cables, use hubs between approved high-speed cables.



- 2. Never attempt to make your own USB cable, or modify one. USB is sensitive to cable impedances and cable lengths. Only use a certified USB cable from a reliable supplier, never a cheap brand or 'something the workshop whipped up'. Your PowerLab is supplied with the proper USB cable. If you need additional USB cables, you should buy high-speed cables (fully shielded, twistedpair, and with standard USB connections) for reliable results.
- 3. Don't disconnect the PowerLab while a program is using it. Disconnecting a PowerLab while it is sampling is a bad idea and may cause problems in some circumstances. It is better to be safe than sorry, and not even disconnect a PowerLab while Chart or Scope is open. Stop sampling and quit the program first before disconnecting a PowerLab from USB.

You can safely turn on or off, or disconnect or reconnect, a USBconnected PowerLab while the computer remains on as long as the application program (Chart or Scope) is off when you do it.

The SCSI Connection

This section is essential reading before trying to use your PowerLab with SCSI. The user's guide for your computer may have a section on connecting SCSI devices, in which case you should read that as well.


The PowerLab is fitted with a SCSI port, which provides an alternative to the USB connection. SCSI (small computer system interface) is simply a communication standard for connecting a single chain of SCSI devices together on a common high-speed bus (information-carrying pathway). A SCSI connection can easily handle the 4 megabytes per second maximum data transfer rate of the PowerLab, although the speed of your computer may limit this. The PowerLab must be placed at the end of the SCSI chain (if it is the only device, then connected directly to the SCSI port of, or to an installed SCSI card in, the computer).

SCSI Connection Rules

 Never alter the SCSI chain while any SCSI devices are on. Disconnecting a SCSI device while there is SCSI hard disk activity

Figure 2–3

A SCSI chain: up to six (or seven) devices can be linked to the computer

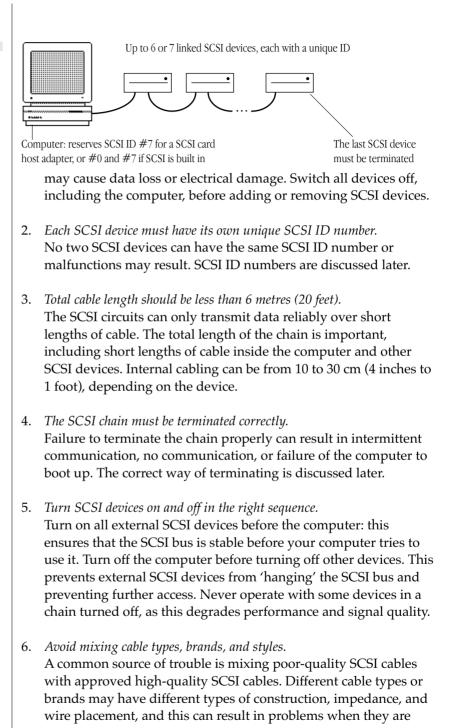

used together. Also, whenever possible, avoid using cable joiners and junctions to extend cable lengths.

7. Never attempt to make your own SCSI cable, or modify one. SCSI is sensitive to cable impedances and cable lengths. Only use an approved SCSI cable from a reliable supplier, never 'something the workshop whipped up'. If you need additional SCSI cables, you should buy cables that are double-shielded (foil and braid), with twisted pairs and shielded connector hoods for the most reliable results. ADInstruments sells approved SCSI cables

Setting the SCSI ID Number

Your PowerLab was set to a SCSI ID number of 4 at the factory before shipping. To change its SCSI ID, turn off the PowerLab, and all other connected devices in the SCSI chain, before changing the SCSI ID number. The SCSI ID switch is located on the back panel of the PowerLab. A small panel with push buttons above and below it displays the SCSI ID number. Press the [–] button above the indicator to decrease it or the [+] button below the indicator to increase it.

Do not set the SCSI ID to numbers already used by other SCSI devices in the chain. The maximum number of devices within the same SCSI chain is eight. Each device in the chain must have a unique SCSI identification number assigned to it. The numbers range from 0 to 7, with 7 the highest priority device. A computer with built-in SCSI usually reserves the values 0 and 7. On the Macintosh, 7 is used for the computer and 0 may be used for an internal SCSI hard disk (an internal SCSI CD, if there is one, usually takes ID 3, as well). On a PC, SCSI ID number 7 is usually reserved by a SCSI card host adapter or built-in SCSI controller, and SCSI ID number 0 is used by an internal SCSI hard disk, if there is one. To prevent most problems, therefore, the numbers 0 or 7 cannot be used for an SP series PowerLab.

Externally connected devices should have readable SCSI ID switches you can check, although some may have ID switches that are difficult to see or read. Internally connected devices should be described in the manual that came with your computer. If you are unsure about the SCSI devices connected to a Macintosh computer, you can use tools such as SCSIProbe (a free utility available from the Internet and user groups), which lists all SCSI devices attached, what they are, and which SCSI ID they use. If you are unsure about the SCSI devices

Figure 2–4 The SCSI ID switch



connected to a Windows computer (particularly internal ones), your PC technical support person or IS manager may be able to help.

To confirm that you have set the right SCSI ID number, turn on the PowerLab and watch to see if the Status indicator flashes yellow a number of times equal to its SCSI ID number.

SCSI Termination Rules


Terminating the SCSI bus is essential to preserve the integrity of the high-speed signals present on the bus. Termination helps keep these signals free from noise and prevents unwanted signal reflections up and down the bus. Without correct termination, the SCSI signals could be swamped by reflections from previous signals, resulting in a jumble of meaningless electronic noise. The way in which you terminate the SCSI chain will depend on the number of SCSI devices you wish to connect to your computer along with the PowerLab. The SCSI connection between your PowerLab and computer is a very good way to provide safe and reliable high-speed data transfer, provided that you follow the rules for proper termination.

A SCSI chain is correctly terminated by placing terminators at each end of the chain. The computer is terminated internally, so forming one end of the chain. (A SCSI host adapter will be terminated: if the SCSI controller and internal SCSI hard disk are both terminated, you may have to alter things — see the documentation that came with the card or computer.) The PowerLab should be at the other end of the chain, with its SCSI termination switch on.

Connecting the PowerLab: Macintosh

Those Macintosh computers without USB (and even some with USB) should have SCSI built in or available through a card. Not all SCSI cards are suitable, especially high-end ones (ultra-wide SCSI and the like). SCSI cards and cables with which the PowerLab should work reliably are sold by ADInstruments. Make sure that you have the correct, 25-pin to 25-pin SCSI cable to connect up your PowerLab. (Most PowerBooks and Duo Docks have a special space-saving square 29-pin SCSI connector, which requires another connector or adapter. This can be purchased separately from your Apple dealer.)

If the PowerLab is the only SCSI device to connect to the computer, then simply connect the two together with the correct SCSI cable, and turn the SCSI termination switch on. If you wish to use multiple SCSI devices with the computer, the first and last devices in the chain should be terminated. The PowerLab must be at the end of the chain, since it has only one SCSI connector, and its internal terminator must be switched on.

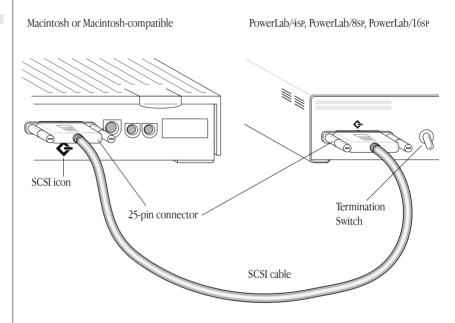

Connecting the PowerLab: Windows


Most modern PC computers should have SCSI available through a card (some have it built in). Not all SCSI cards are suitable, especially high-end ones (ultra-wide SCSI and the like). ADInstruments sells SCSI cards and cables with which the PowerLab should work reliably. (It may not be possible to find a cable to fit all connections, though.)

Make sure that you have the correct SCSI cable to connect up your PowerLab. It requires a 25-pin D-style connector for the PowerLab. If you have a portable computer with a PC card SCSI adapter, then the card probably needs the specific cable that came with that particular card. You may have to purchase separately a cable adapter to connect the end of your SCSI cable to the 25-pin connector for the PowerLab.

Figure 2–5

Connecting a PowerLab to the computer (look for the SCSI icon if the back panel is different) The card should have full instructions on installation with it. The SCSI adapter card will usually be terminated, and at one end of the chain (this is always the case for PC card SCSI adapters), although if the computer has built-in SCSI, you may have an internal SCSI device in your PC that is terminated, with the host adapter in the middle of the chain. If there is a problem with cables, or you are unsure about any devices, such as a hard disk, connected to your computer on SCSI internally, your PC technical support person or IS manager should be able to help you.

If the PowerLab is the only SCSI device to connect to the computer, then simply connect the two together with the correct SCSI cable, and turn the SCSI termination switch on. If you wish to use multiple SCSI devices with the computer, the first and last devices in the chain should be terminated. The PowerLab must be at the end of the chain, since it has only one SCSI connector, and its internal terminator must be switched on.

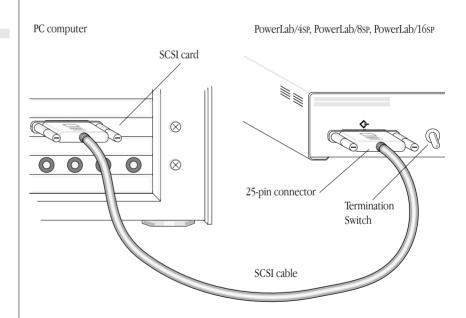

SCSI Cards with Windows NT or Older Computers

Windows 95 and 98 both come with a number of generic SCSI drivers already installed that should work with most cards, including any that ADInstruments supplies. Windows NT may not have the right

Figure 2–6 Connecting a PowerLab to the computer (25-pin connector shown) SCSI driver available, so the driver might have to be installed manually, perhaps from the original NT system CD. The SCSI card should have instructions on the driver it needs and how to install it.

Older, pre-PCI computers require a special ISA or ISO bus SCSI card. Such a card should come with its own manual, and may take some effort to set up. It will probably require assistance from a PC technical support person.

SCSI Strikes Back

Not following the SCSI device connection and termination protocols can lead to more problems than you might think. The possible dangers include loss of data, intermittent communications, corruption of data on a SCSI hard disk, failure of the computer to boot from a SCSI drive, and failure of the computer to recognise SCSI devices connected to it. Such problems are very unlikely, however, if you follow the rules.

A P P E N D I X A

A

Technical Aspects

This appendix describes some of the important technical aspects of the PowerLab/4SP, PowerLab/8SP, and PowerLab/16SP, to give some insight into how they work. You do not need to know the material here to use your PowerLab. It is likely to be of especial interest to the technically minded, indicating what the PowerLab can and cannot do, and its suitability for particular purposes. You should not use it as a service manual: remember that user modification of the PowerLab voids your rights under warranty.

How Does it Work?

The PowerLab is essentially a smart peripheral device specifically designed to perform the various functions needed for data acquisition, signal conditioning, and pre-processing. It contains its own microprocessor, memory, and specialised analog amplifiers for signal conditioning. The block diagrams in Figure A–1, Figure A–2, and Figure A–3 show the essential elements of each PowerLab.

All sampling, output, and communication functions are controlled by an internal 32-bit, 68340 microprocessor running at 16 MHz (16 million cycles per second). This CPU (central processing unit) has access to 512 kilobytes of RAM (random access memory) for data storage and buffering prior to transmission to the computer. The PowerLab makes use of FIFO (first in, first out) memory devices, which provide a temporary storage area for both analog input and analog output signals. These storage areas help increase the processing capacity of the hardware by relieving the load on the PowerLab microprocessor. The PowerLab can use SCSI or USB, whether built-in or via a card, to communicate with the computer. SCSI (small computer system interface) provides data transfer rates of up to 4 MB per second on fast computers. The lower-speed USB (universal serial bus) provides data transfer rates of up to 500 kB per second.

The PowerLab has four (PowerLab/4SP), eight (PowerLab/8SP), or 16 (PowerLab/16SP) analog inputs, used to record external signals. The outputs of their input amplifiers are multiplexed to a 16-bit ADC


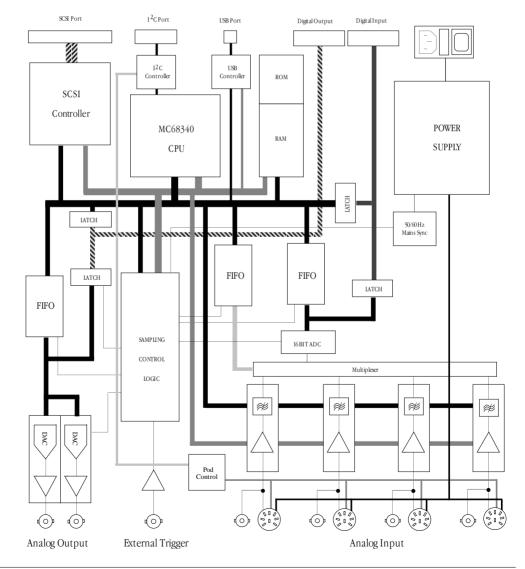


Figure A–1 Block diagram of the PowerLab/4SP (analog-to-digital converter). The ADC can sample at up to 200,000 samples per second. The CPU assembles groups of samples into blocks and then transmits them to the computer, where the application program receives, records, and displays the data. Sampled data from the ADC are stored in FIFOs until the CPU is ready to read them. The FIFOs let the CPU know when they are getting full, at which point it will read all of their data very quickly. This frees the CPU from the need to read every sample from the ADC as it is ready, so that it can perform other tasks.


The external trigger input (marked 'Trigger' on the front panel) is connected to a comparator circuit that triggers when the input



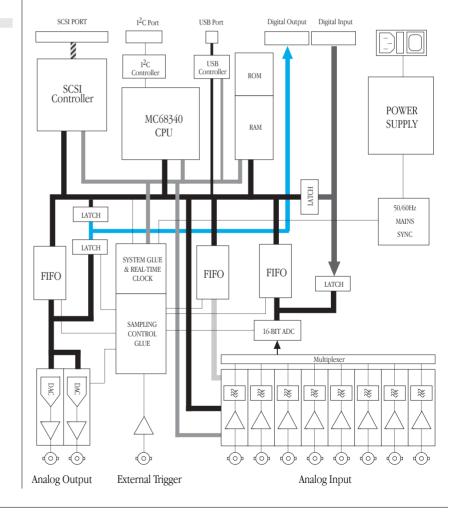

voltage exceeds 2.9 volts. This signal is fed to circuitry that notifies the CPU that an external trigger event has been detected. The CPU then carries out the task for which the trigger is being used (such as pre-triggering or post-triggering). When the trigger threshold is crossed, the indicator beside the trigger connector glows yellow.

Two 14-bit DACs (digital-to-analog converters) are used to control the analog outputs of the PowerLab (marked 'Output 1' and 'Output 2' on the front panel). The DACs produce waveforms under software control that are fed through an attenuation network to produce different full-scale ranges. The signals are then buffered by a power

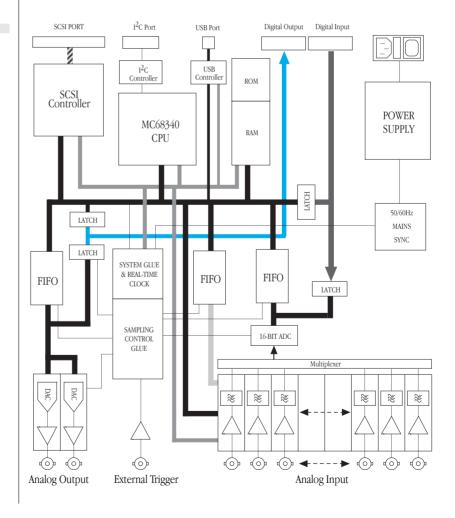



▲ Caution

PowerLab inputs and outputs are not electrically isolated. Human subjects must not be connected directly to the PowerLab. If such measurements are to be made, an isolated front-end must be used, such as the Bio Amp. amplifier to give them the capability to drive heavier loads. The PowerLab can store stimulation data in FIFOs, so that it can stimulate at a frequency independent of the analog input sampling rate. The FIFOs also allow the generation of much faster stimuli and continuous waveforms, with little or no load on the CPU.

The PowerLab is also fitted with an I^2C front-end expansion port. This 9-pin port supplies both power and control to ADInstruments front-ends using a 4-wire serial bus (two wires for standard I^2C and two control lines).

The digital input and output connectors provide a means of using the PowerLab to monitor and to control external devices respectively. The digital input monitors state changes: the eight lines of the connector allow monitoring of up to eight devices. The digital output can turn on and off external devices or can signal to some other device. The eight lines of the connector allow control of up to eight devices. These recording automation and control features are available in the Macintosh version of Chart, but not in the current version of Chart for Windows (v3.4), or in Scope.

The PowerLab uses an IEC601-1 compliant linear power supply, not a switching power supply, which makes it a little heavier, but a lot quieter electrically. It is important to note that the PowerLab has a limited amount of power available for external devices. Because of these power restrictions, you should not use the PowerLab as a power source for external devices other than front-ends produced by ADInstruments.

The Analog Inputs


PowerLab input amplifiers have been designed with a considerable amount of computer-controlled gain (up to × 5000). Thus it is possible to record a variety of signals without any external pre-amplification. Each analog input is a separate DC amplifier, with programmable gain able to be set independently (the gain is set through the software range control: the less the range, the more the gain). The input amplifiers of the PowerLab/8SP and PowerLab/16SP are singleended, but can be set by the software to be either positive or negative (inverting or non-inverting). The PowerLab/4SP input amplifiers can be set by the software to be either single-ended — positive or negative — or differential. In the differential setting, the amplifier

▲ Caution

Applying more than ± 15 V to the input can damage the channel input circuits.

measures the difference between the positive and negative inputs of a Pod connector, irrespective of ground. Differential signals can only be recorded using the Pod connectors on the PowerLab/4sP; the BNC analog inputs on the PowerLabs are all single-ended.

It is important to note that the PowerLab grounds the inputs to amplifiers not in use. It also grounds each amplifier and measures the DC offset voltage when the gain is changed. In this way, the software corrects for any DC drift or offset in the circuits that may develop over time or between readings. The operation of the input amplifiers is illustrated by the block diagrams below.

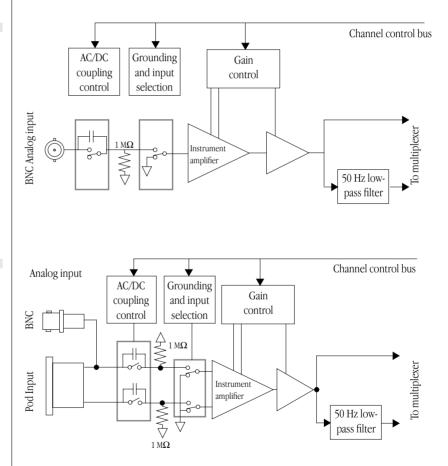

The input amplifiers can be set to pass both DC and AC signals, or to pass only AC signals without passing DC signals. Input impedance is one megohm (1 $M\Omega$).

Figure A-4

Block/schematic diagram of the input amplifier circuitry for BNC-only inputs

Figure A-5

Block/schematic diagram of the input amplifier circuitry for dual BNC and Pod inputs

Each input produces either a filtered or an unfiltered signal. Each input is fitted with a software-selectable low-pass filter (a fourthorder anti-aliasing filter), which can be set at a range of frequencies from 20 kHz down to 1 kHz. The unfiltered signal has an effective bandwidth of 45 kHz (within the Nyquist frequency when sampling at 100,000 samples per second).

PowerLab Accuracy

Your PowerLab was calibrated at the factory to an accuracy of better than 0.1%. Some 'zero drift' or 'gain drift' can occur with time. This can affect the accuracy of measurements, especially at the highest input gains. The unit can be recalibrated, but in most circumstances this is not necessary in its lifetime. There are several reasons for this.


DC drift compensation. Each time that recording is started manually or by triggering or the gain is changed (that is, very often in most cases), the input to the amplifier is grounded and any DC due to the amplifier's drift with temperature and age is measured. The measured voltage is removed from the readings for that input through software correction, in a process transparent to the user.

Calibration facilities. It is recommended and sound practice to calibrate a measuring system from the transducer to the output. After applying two known values to a transducer (say at 20% and 80% of full scale) and recording the signal, you can use the units conversion feature of ADInstruments applications to convert and display transducer readings in the appropriate units. This will compensate for any minor inaccuracies in amplifier gain and transducer calibration.

The External Trigger


The external trigger provides a digital input for synchronising sampling to external devices. The external trigger input represents 1 TTL (transistor–transistor logic) load. The input is off for input voltages between –12 V and the external trigger level, and on between that and +12 V. The input will be overloaded if the voltage is outside the range –12 V to +12 V. The external trigger level is 2.9 V \pm 0.25V; once on, the trigger turns off at 1.8 V \pm 0.25V, giving a hysteresis voltage of 1.1 V. When the external trigger input is on, the indicator beside it is on.


The PowerLab has an optically isolated external trigger input, thus there is no direct connection between the external trigger ground and the ground of the device connected to it. This removes ground noise and current problems, and improves static discharge immunity. In order for the external trigger to work, though, a voltage must be applied between the outer ring and the inner pin on the connector. Applying a voltage to the centre pin won't result in triggering (as it could on earlier PowerLabs). The equivalent circuit of the external trigger is basically an LED (light-emitting diode) in series with a resistor, thus the device driving the external trigger must be capable of supplying at least 1 mA of current to cause the optocoupler to work. (The second diode shown below protects the optocoupler.)

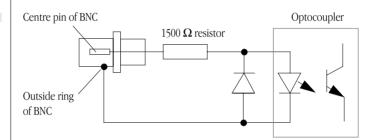

The Analog Output

The analog outputs provide computer-controlled variable outputs (±10 V) that can be used with the Chart and Scope applications either directly as a stimulator, or to control peripheral devices. All stimulation voltage is generated by the PowerLab via the output sockets on the front of the PowerLab (marked Output 1 and Output 2), giving positive, negative, differential, or independent stimuli, depending on the sockets used and the software settings.

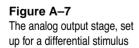

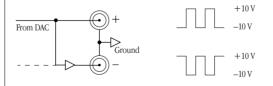

By default, the outputs are used for complementary (differential) stimulation, where Output 1 is positive and Output 2 is negative. When Output 1 is used, a positive stimulus voltage (set up in an application) gives a positive voltage output, and a negative voltage a

Figure A–6 The equivalent circuit of the external trigger input

negative one. When Output 2 is used, the voltage outputs are inverted. When both output sockets are used, the stimulus is the difference between the voltages at the positive and negative outputs: you could generate up to a 20-volt pulse, given a ± 10 V stimulus.

The outputs can be used for independent stimulation in Scope and the Macintosh version of Chart, but not in the current version of Chart for Windows (v3.4). The details are discussed in the software user's guides.

Connections

This section of the appendix contains 'pinout' and electrical details of some of the connectors fitted to the PowerLab. You should read it carefully before attempting to connect cables other than those supplied with the unit to the PowerLab. Using cables that are wired incorrectly can cause internal damage to the PowerLab and will void your rights under warranty.

USB Port

The PowerLab is fitted with a USB (universal serial bus) port, which provides an alternative to the SCSI connection. This works with computers with USB connectors or a USB card. USB needs a Power Macintosh with Mac OS 8.5 or later, or a PC with Windows 98 or later.



Your PowerLab is supplied with a proper high-speed USB cable. If you need additional USB cables, you should buy high-speed cables (fully shielded, twisted-pair, and with standard USB connections) for reliable results.


SCSI Port

The PowerLab uses a single 25-pin D-type female connector for the SCSI port. If you are using a Macintosh, you probably need a 25-pin

Figure A–8 The USB connector

to 25-pin cable. If you are using a Windows PC, the connection at the other end of the cable depends on your set-up.

Figure A–9 The 25-pin SCSI connector

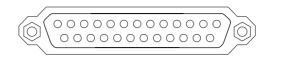

SCSI is sensitive to cable impedances and connection lengths. The incorrect wiring of a SCSI cable can result in damage to the SCSI circuitry, or random SCSI activities being performed. You should, therefore, use only approved SCSI cables such as the ones available from ADInstruments, and never attempt to make your own.

I²C Expansion Port

The I²C port on the back panel of the PowerLab provides expansion support for ADInstruments front-ends. This port provides both power and control signals for these front-ends. The I²C bus has a daisy-chain structure that allows simple connection of additional front-ends to the system. A PowerLab can have as many front-ends connected to it as it has analog inputs, roughly speaking. You should not attempt to run other external devices from the I²C port: it is designed for use only with ADInstruments front-ends. Only 50 mA maximum current can be provided through this bus, so it should not be used for third-party devices drawing more current.

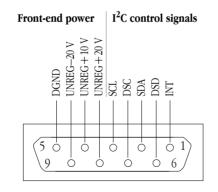



Figure A–10 The pin assignments for the I^2C port G

Digital Input and Output Ports

The digital input port and digital output port are 15-pin connectors situated on the back panel of the PowerLab. The eight digital input lines respond to standard TTL signals with a threshold of 2.2 V. The eight digital output lines can turn on and off, or signal to, up to eight external TTL devices. The digital output lines are capable of driving five TTL loads overall (50 mA maximum load per line; 250 mA maximum total load).

Digital input



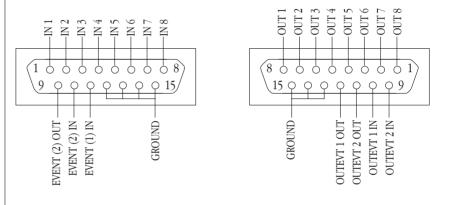

These recording automation and control features are available in the Macintosh version of Chart, but not in the current version of Chart for Windows (v3.4), or in Scope.


Pod Connectors


The Pod connectors on the PowerLab/4SP are 8-pin DIN connectors. (The PowerLab/8SP and PowerLab/16SP have single-ended BNC inputs only.) Do not attempt to record from both the BNC and Pod connectors for an input at the same time, or the signals will compete.

Pod connectors allow the connection of ADInstruments Pods small, low-cost units that provide alternatives to front-ends for specific tasks, for use with precalibrated transducers and so on. (Pods are supported by Scope v3.6.3 and later, Chart for Macintosh v3.6.3 and later, and Chart for Windows v3.4.7 and later.) The Pod

Figure A–11 The pin assignments for the digital input and output connectors

connectors on the PowerLab/4SP do not handle transducers directly unless the transducers are so labelled (unsuitable transducers will give a very weak signal). Transducers designed for direct connection can be provided with power and control, since the Pod connectors provide some functions of the I²C output as well as alternative analog inputs to the BNC connectors.

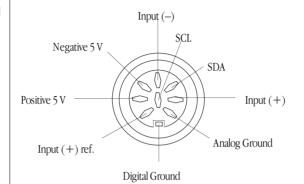

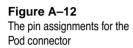

A P P E N D I X B

B

Troubleshooting

This appendix describes most of the common problems that can occur with your PowerLab recording unit. It covers how these problems are caused, and what you can do to alleviate them. If the solutions here do not work, your software guide may contain remedies for possible software problems. If none of the solutions here or in the software guide appears to help, then consult your ADInstruments distributor.

Problems: Macintosh

Nearly all the problems that users encounter are connection problems and SCSI problems. Improper connections can result in a loss of all or some of a signal. SCSI difficulties can cause data loss, failure of the computer to recognise that SCSI devices are attached to it, and difficulties using other SCSI devices. Incompatible SCSI adapter cards and cables may also cause problems. Obviously, the SCSI problems will not apply if you use USB.

The PowerLab Status indicator light flashes yellow then red when the PowerLab is turned on

If the pattern is simply an alternating red flash and yellow flash, then the SCSI ID has been set to the numbers 0 or 7, which cannot be used by an SP series PowerLab. Only 1 to 6 are valid SCSI ID numbers.

• Switch off your PowerLab and any other connected devices, change the SCSI ID for the PowerLab using the control on the back panel, wait about five seconds, and then switch it back on. Check any other connected SCSI devices as well at the same time.

▼ **Refer** Self-Test, p. 16 An internal problem has been discovered by the diagnostic self-test (this is performed by the PowerLab each time it powers up).

• Turn everything off, and then after at least five seconds turn the PowerLab back on again. This should clear a temporary problem. If not, then the PowerLab may need repair. Take note of the flashing pattern, and consult your ADInstruments distributor.

The computer refuses to boot with the PowerLab connected, or the computer can't find the PowerLab

The PowerLab is off or the power is switched off at the wall, the power cable is not connected firmly, or a fuse has blown.

• Check switches, power connections, and fuses.

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

The SCSI adapter card is incompatible with the computer or the PowerLab or both. We cannot guarantee that the PowerLab will work reliably with a card we did not supply.

• Try another card, preferably one we supplied.

The PowerLab has an internal problem or has 'hung'.

• Turn everything off, and then after at least five seconds turn the PowerLab back on again. Turn on the computer and try using the software again.

The PowerLab is connected via USB and the computer does not properly support USB.

 USB needs a Power Macintosh with Mac OS 8.5 or later; it simply will not work with earlier hardware, such as a 68K Macintosh, or earlier operating systems, such as System 7: use SCSI with these.

The computer hangs up while recording, or there is data loss

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

The PowerLab doesn't work or the program crashes after a short time

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

The SCSI adapter card is incompatible with the computer or the PowerLab or both. We cannot guarantee that the PowerLab will work reliably with a card we did not supply.

• Try another card, preferably one we supplied.

Partitioning problems on the hard disk (while rare, this can happen even with new disks).

• Use disk-checking utility applications to check the SCSI driver and root directory on the hard disk. Partitioning programs may

also have effective low-level disk checking. The programs will probably fix the problems, or otherwise indicate what needs to be done (such as reinstalling the SCSI driver), and their manuals should explain the technical details involved.

The PowerBook computer stops working normally, and instead shows a diamond with its SCSI ID number on screen

The PowerBook is acting as a SCSI hard disk. This means that the wrong (30-pin) connector and cable were used to connect it to the PowerLab. You need a different (29-pin) cable and connector to make a normal SCSI connection to peripherals.

• Use the correct cable to connect your PowerBook to the PowerLab (the cable should be available from your Apple dealer).

Signals seem weak or interacting on the PowerLab/4SP

You may be using both the BNC connector and the Pod connector for an input at the same time, with resultant signal competition.

• Make certain that you use only *one* of the alternative connectors (BNC or Pod) for an input when recording.

Problems: Windows

Most problems that users encounter are connection problems and SCSI problems. Improper connections can result in a loss of all or some of a signal. SCSI difficulties can cause data loss, failure of the computer to recognise that SCSI devices are attached to it, and difficulties using other SCSI devices. Incompatible SCSI adapter cards and cables may also cause problems. Obviously, the SCSI problems will not apply if you use USB.

The PowerLab Status indicator light flashes yellow then red when the PowerLab is turned on

If the pattern is simply an alternating red flash and yellow flash, then the SCSI ID has been set to the numbers 0 or 7, which cannot be used by an SP series PowerLab. Only 1 to 6 are valid SCSI ID numbers. • Switch off your PowerLab and any other connected devices, change the SCSI ID for the PowerLab using the control on the back panel, wait about five seconds, and then switch it back on. Check any other connected SCSI devices as well at the same time.

An internal problem has been discovered by the diagnostic self-test (this is performed by the PowerLab each time it powers up).

• Turn everything off, and then after at least five seconds turn the PowerLab back on again. This should clear a temporary problem. If not, then the PowerLab may need repair. Take note of the flashing pattern, and consult your ADInstruments distributor.

When Windows starts up, it doesn't recognise the PowerLab

This should happen only the first time PowerLab hardware is connected to the computer. Windows may bring up the New Hardware wizard, and ask if you want to install a driver.

• Leave the wizard on its default settings and click the OK button. Insert the CD, or Disk 1 of the provided installation disks. The PowerLab setup information file should be selected in the dialog box that appears. Click the OK button, and continue.

There is an alert about ASPI when the software starts up

ADInstruments software needs ASPI (advanced SCSI programming interface) to communicate with the PowerLab. An alert appears if it can't be found. ASPI is present by default in Windows 95 or later (although it could have been removed), but not in Windows NT 4. A version suitable for most SCSI cards supplied by ADInstruments is provided on the software CD. Other SCSI cards should come with instructions on what to do.

• Install the appropriate version of ASPI on the computer, or ask your PC technical support person or IS manager to do it.

The computer refuses to boot with the PowerLab connected

The SCSI adapter card is incompatible with the computer or the PowerLab or both. We cannot guarantee that the PowerLab will work

▼ **Refer** Self-Test, p. 16 reliably with a card we did not supply, and some computers and SCSI adapter cards just won't work together.

• Try another card, preferably one we supplied. Ask your PC technical support person or IS manager to check compatibility.

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

The computer is running off a SCSI hard drive, and the PowerLab has an internal problem or has 'hung'.

• Turn everything off, and then after at least five seconds turn the PowerLab back on again. Turn on the computer and try using the software again.

The computer can't find the PowerLab

The PowerLab is off or the power is switched off at the wall, the power cable is not connected firmly, or a fuse has blown.

• Check switches, power connections, and fuses.

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

The PowerLab has an internal problem or has 'hung'.

• Turn everything off, and then turn the PowerLab back on again after at least five seconds. Turn on the computer and try using the software again.

Windows NT may not have the right SCSI driver available (it does not have many already installed, as do Windows 95 and 98).

• The driver might have to be installed manually, perhaps from the original NT system CD. The SCSI card should have instructions on the driver it needs and how to perform the installation.

The PowerLab is connected via USB and the computer does not properly support USB.

• USB needs a USB 1.1-compliant PC with Windows 98, Me, 2000 or later; it simply will not work with earlier operating systems, such as Windows 95 or Windows NT 4: use SCSI with these.

The computer hangs up while recording, or there is data loss

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

The PowerLab doesn't work or the program crashes after a short time

The SCSI adapter card is incompatible with the computer or the PowerLab or both. We cannot guarantee that the PowerLab will work reliably with a card we did not supply, and some computers and SCSI adapter cards just won't work together.

• Try another card, preferably one we supplied. Ask your PC technical support person or IS manager to check compatibility.

A poor connection between PowerLab and computer, or bad cable.

• Ensure that the cable is firmly attached at both ends and try again. If there is still a problem, try a new cable.

SCSI problems: two SCSI devices on the bus have the same SCSI ID, or termination is incorrect or unsuitable.

• Check that each device has a unique ID number and that the SCSI chain is terminated correctly. (As a last resort, if the PowerLab is the only connected SCSI device, try leaving it unterminated.)

Signals seem weak or interacting on the PowerLab/4SP

You may be using both the BNC connector and the Pod connector for an input at the same time, with resultant signal competition.

• Make certain that you use only *one* of the alternative connectors (BNC or Pod) for an input when recording.

A P P E N D I X C

Specifications

PowerLab/4sp Specifications

Input

C

Number of inputs:	4
Input configuration:	Single-ended or differential (the latter only through the Pod connectors)
Amplification ranges:	$\pm 2 \text{ mV}$ to $\pm 10 \text{ V}$ full scale in 12 steps
	$\begin{array}{c} \pm 10 \text{ V} \\ \pm 5 \text{ V} \\ \pm 2 \text{ V} \\ \pm 1 \text{ V} \\ \pm 0.5 \text{ V} \\ \pm 0.2 \text{ V} \\ \pm 0.1 \text{ V} \\ \pm 50 \text{ mV} \\ \pm 20 \text{ mV} \\ \pm 10 \text{ mV} \\ \pm 5 \text{ mV} \\ \pm 2 \text{ mV} \end{array}$
Maximum input voltage:	±15 volts
Input impedance:	$\approx 1 \text{ M}\Omega \parallel 47 \text{ pF } @ \text{ DC}$
Low-pass filtering:	1, 2, 5, 10, or 20 kHz (software-selectable)
AC coupling:	DC or 0.1 Hz (software-selectable)
Frequency response (-3 dB):	20 kHz @ ±10 V full scale, all ranges

DC drift:	Software-corrected
CMRR (differential):	96 dB @ 50 Hz (typical)
Input crosstalk:	–110 dB typical
Input noise:	$<1\mu V_{rms}$ referred to input
Pod connectors:	Combine power, I ² C and single-ended or differential analog input signals on one connector, support Pods, particular transducers, etc.
Supply voltage: Maximum current: Communications: Signal input: Connector type:	±5 V regulated 50 mA per pod port 2-wire I ² C Positive and negative analog inputs 8-pin DIN.
Sampling	
ADC resolution:	16 bits (150 μV resolution at 10 V range)
Linearity error:	±1 LSB (from 0 °C to 70 °C)
Maximum sampling rates:	200 kHz on one channel; 100 kHz on two channels; 40 kHz on three or four channels.
Available sampling rates:	200 kHz down to 0.003 Hz using Chart 200 kHz down to 2 Hz using Scope 200 kHz to 1 Hz using Chart for Windows
	Sampling on USB rather than SCSI may limit the maximum sampling rates.
Output Amplifier	
Number of outputs:	2
Output configuration:	Single-ended or differential
Output resolution:	14 bits (0.6 mV resolution at 10 V range)
Maximum output current:	±25 mA
Output impedance:	0.1Ω typical
Slew rate:	10 V/µs
Settling time:	$2\mu s$ (to 0.01% of FSR for LSB change)

Linearity error:	±1 LSB (from 0 °C to 70 °C)	
Output ranges:	$\pm 200~\text{mV}$ to $\pm 10~\text{V}$ full scale in six steps	
	± 10 V ± 5 V ± 2 V ± 1 V ± 500 mV ± 200 mV	
External Trigger		
Trigger threshold:	+2.9 V (TTL compatible)	
Hysteresis:	1.1 V (turns off at +1.8 V)	
Input load:	1 TTL load	
Maximum input voltage:	±12 V	
Minimum pulse width:	5 µs	
Microprocessor and Data Communication		
Processor:	16 MHz 68340	
Memory:	512 K RAM	
Data communication:	SCSI (up to 4 megabytes/s, maximum, dependent on computer)	

Expansion Ports

I ² C expansion port:	Power and control bus for front-end units. Supports up to 16 front-ends, but limited to the PowerLab's free connectors. Interface communications rate of up to 10,000 bits/s.
Digital output:	8 independent lines, TTL output level (50 mÅ maximum load per line; 250 mA maximum total load)
Digital input:	8 independent lines, TTL input level, threshold 2.2 V

USB (up to 500 kilobytes/s, maximum dependent on setup)

Physical Configuration

Dimensions (w \times h \times d):	300 mm × 60 mm × 300 mm (11.8" × 2.4" × 11.8")
Weight:	4.5 kg (9 lb 13 oz)
Operating voltage:	220-240 V or 110-120 V (internally set)
Nominal power needs:	12.5 VA (no front-ends attached)
Maximum power needs:	18 VA (full complement of front-ends)
Replacement fuses:	100–120 V: 1 Amp (5 × 20 mm) slow blow (voltage rating 120 or 250 V)
	220–240 V: 400 mA (5 × 20 mm) slow blow (voltage rating 250 V only)
Operating temperature range:	0 to 35 °C, 0 to 90% humidity (non- condensing)

PowerLab/8sp Specifications

Input

Number of inputs:	8
Input configuration:	Single-ended
Amplification ranges:	$\pm 2~mV$ to $\pm 10~V$ full scale in 12 steps
	$\begin{array}{c} \pm 10 \text{ V} \\ \pm 5 \text{ V} \\ \pm 2 \text{ V} \\ \pm 1 \text{ V} \\ \pm 0.5 \text{ V} \\ \pm 0.2 \text{ V} \\ \pm 0.1 \text{ V} \\ \pm 50 \text{ mV} \\ \pm 20 \text{ mV} \\ \pm 10 \text{ mV} \\ \pm 5 \text{ mV} \\ \pm 2 \text{ mV} \end{array}$
Maximum input voltage:	±15 volts
Input impedance:	$\approx 1 \ M\Omega \parallel 47 \ pF \ @ \ DC$

Low-pass filtering:	1, 2, 5, 10, or 20 kHz (software-selectable)
AC coupling:	DC or 0.1 Hz (software-selectable)
Frequency response (-3 dB):	20 kHz @ ±10 V full scale, all gain ranges
DC drift:	Software-corrected
CMRR (differential):	96 dB @ 50 Hz (typical)
Input crosstalk:	–110 dB typical
Input noise:	$<1\mu V_{rms}$ referred to input
Sampling	
ADC resolution:	16 bits (150 μV resolution at 10 V range)
Linearity error:	±1 LSB (from 0 °C to 70 °C)
Maximum sampling rates:	200 kHz on one channel; 100 kHz on two channels; 40 kHz on three or four channels; 20 kHz on five to eight channels.
Available sampling rates:	200 kHz down to 0.003 Hz using Chart 200 kHz down to 2 Hz using Scope 200 kHz to 1 Hz using Chart for Windows
	Sampling on USB rather than SCSI may limit the maximum sampling rates.
Output Amplifier	
Number of outputs:	2
Output configuration:	Single-ended or differential
Output resolution:	14 bits (0.6 mV resolution at 10 V range)
Maximum output current:	±25 mA
Output impedance:	0.1Ω typical
Slew rate:	10 V/µs
Settling time:	$2\mu s$ (to 0.01% of FSR for LSB change)
Linearity error:	±1 LSB (from 0 °C to 70 °C)

Output ranges:	$\pm 200 \text{ mV}$ to $\pm 10 \text{ V}$ full scale in six steps
	± 10 V ± 5 V
	± 2 V ± 1 V
	$\pm 500 \text{ mV}$ $\pm 200 \text{ mV}$
External Trigger	
Trigger threshold:	+2.9 V (TTL compatible)
Hysteresis:	1.1 V (turns off at +1.8 V)
Input load:	1 TTL load
Maximum input voltage:	±12 V
Minimum pulse width:	5 μs
Microprocessor and Dat	a Communication
Processor:	16 MHz 68340
Memory:	512 K RAM
Data communication:	SCSI (up to 4 megabytes/s, maximum, dependent on computer)
	USB (up to 500 kilobytes/s, maximum dependent on setup)
Expansion Ports	
I ² C expansion port:	Power and control bus for front-end units. Supports up to 16 front-ends, but limited to the PowerLab's free connectors. Interface communications rate of up to 10,000 bits/s.
Digital output:	8 independent lines, TTL output level (50 mÅ maximum load per line; 250 mA maximum total load)
Digital input:	8 independent lines, TTL input level, threshold 2.2 V

Physical Configuration

Dimensions (w \times h \times d):	300 mm × 60 mm × 300 mm (11.8" × 2.4" × 11.8")
Weight:	4.8 kg (10 lb 8 oz)
Operating voltage:	220–240 V or 110–120 V (internally set)
Nominal power needs:	17 VA (no front-ends attached)
Maximum power needs:	30 VA (full complement of front-ends)
Replacement fuses:	100–120 V: 1 Amp (5 × 20 mm) slow blow (voltage rating 120 or 250 V)
	220–240 V: 400 mA (5 \times 20 mm) slow blow (voltage rating 250 V only)
Operating temperature range:	0 to 35 °C, 0 to 90% humidity (non- condensing)

PowerLab/16sp Specifications

Input

Number of inputs:	16
Input configuration:	Single-ended
Amplification ranges:	$\pm 2 \text{ mV}$ to $\pm 10 \text{ V}$ full scale in 12 steps
	$\begin{array}{c} \pm 10 \text{ V} \\ \pm 5 \text{ V} \\ \pm 2 \text{ V} \\ \pm 1 \text{ V} \\ \pm 0.5 \text{ V} \\ \pm 0.2 \text{ V} \\ \pm 0.1 \text{ V} \\ \pm 50 \text{ mV} \\ \pm 20 \text{ mV} \\ \pm 10 \text{ mV} \\ \pm 5 \text{ mV} \\ \pm 2 \text{ mV} \end{array}$
Maximum input voltage:	±15 volts
Input impedance:	$\approx 1 \ M\Omega \parallel 47 \ pF \ @ \ DC$


X (11. 1	
Low-pass filtering:	1, 2, 5, 10, or 20 kHz (software-selectable)
AC coupling:	DC or 0.1 Hz (software-selectable)
Frequency response (-3 dB):	20 kHz @ ±10 V full scale, all gain ranges
DC drift:	Software-corrected
CMRR (differential):	96 dB @ 50 Hz (typical)
Input crosstalk:	–110 dB typical
Input noise:	$< 1 \mu V_{rms}$ referred to input
Sampling	
ADC resolution:	16 bits (150 μV resolution at 10 V range)
Linearity error:	±1 LSB (from 0 °C to 70 °C)
Maximum sampling rates:	200 kHz on one channel; 100 kHz on two channels; 40 kHz on three or four channels; 20 kHz on five to eight channels; 10 kHz on nine to sixteen channels.
Available sampling rates:	200 kHz down to 0.003 Hz using Chart 200 kHz down to 2 Hz using Scope 200 kHz to 1 Hz using Chart for Windows
	Sampling on USB rather than SCSI may limit the maximum sampling rates.
Output Amplifier	
Number of outputs:	2
Output configuration:	Single-ended or differential
Output resolution:	14 bits (0.6 mV resolution at 10 V range)
Maximum output current:	±25 mA
Output impedance:	0.1Ω typical
Slew rate:	10 V/µs
Settling time:	$2\mu s$ (to 0.01% of FSR for LSB change)
Linearity error:	± 1 LSB (from 0 °C to 70 °C)

Output ranges:	$\pm 200 \text{ mV}$ to $\pm 10 \text{ V}$ full scale in six steps
	$\pm 10 V$ $\pm 5 V$ $\pm 2 V$ $\pm 1 V$ $\pm 500 mV$ $\pm 200 mV$
External Trigger	
Trigger threshold:	+2.9 V (TTL compatible)
Hysteresis:	1.1 V (turns off at +1.8 V)
Input load:	1 TTL load
Maximum input voltage:	±12 V
Minimum pulse width:	5 µs
Microprocessor and Da	ta Communication
Processor:	16 MHz 68340
Memory:	512 K RAM
Data communication:	SCSI (up to 4 megabytes/s, maximum, dependent on computer)
	USB (up to 500 kilobytes/s, maximum dependent on setup)
Expansion Ports	
I ² C expansion port:	Power and control bus for front-end units. Supports up to 16 front-ends, but limited to the PowerLab's free connectors. Interface communications rate of up to 10,000 bits/s.
Digital output:	8 independent lines, TTL output level (50 mÅ maximum load per line; 250 mÅ maximum total load)
Digital input:	8 independent lines, TTL input level, threshold 2.2 V

Physical Configuration

z set)
nds)
w blow
ow blow
1-

ADInstruments reserves the right to alter these specifications at any time.

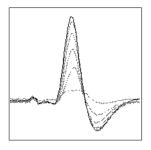

Glossary

This covers terms used in this owner's guide, those for ADInstruments front-ends, and the user's guides for ADInstruments software. General and specific computer terminology should be covered in the material that came with your computer, or in a recent dictionary.

AC coupling. A filter option. When AC coupling is chosen, a 0.1 Hz high-pass filter before the first amplification stage removes DC and frequency components below 0.1 Hz. This removes slowly changing baselines.

ADC (analog-to-digital converter). A device that converts analog information into some corresponding digital voltage or current.

amplitude. The maximum vertical distance of a periodic wave from the zero or mean position about which the wave oscillates.

analog. Varying smoothly and continuously over a range. An analog signal varies continuously over time, rather than changing in discrete steps.

analog input. This refers to the connectors on the front of the PowerLab marked 'Input'. These inputs are designed to accept up to ± 10

volts. Inputs can be either single-sided or differential (the latter only in the case of the Pod connectors).

analog output. This refers to the connectors on the front of the PowerLab marked 'Output'. The analog output provides a software-controlled variable output (±10 V) that can be used with applications either directly as a stimulator, or to control peripheral devices.

analysis. When the PowerLab is not physically connected to the computer, then ADInstruments software can be used to analyse and manipulate existing files if the analysis option is chosen.

BNC. A sort of cable or connector; a BNC-to-BNC cable connects two BNC connectors.

bridge transducer. A type of transducer using a Wheatstone bridge circuit. In its basic form, the bridge consists of four two-terminal elements (usually strain gauges) connected to form a quadrilateral. An excitation source is connected across one diagonal, and the transducer output is taken across the other. bus. A data-carrying electrical pathway.

Chart. An application supplied with a PowerLab that emulates a multi-channel chart recorder, with other powerful options. (Macintosh and Windows versions differ.)

connector. A plug, socket, jack, or port used to connect one electronic device to another (via a cable): a PowerLab to a computer, say.

CPU (central processing unit). A hardware device that performs logical and arithmetical operations on data as specified in the instructions: the heart of most computers.

DAC (digital-to-analog converter). A device that converts digital information into some corresponding analog voltage or current.

DC offset. The amount of DC (direct current) voltage present at the output of an amplifier when zero voltage is applied to the input; or the amount of DC voltage present in a transducer in its equilibrium state.

differential input. Input using both positive and negative inputs on a PowerLab. The recorded signal is the difference between the positive and negative input voltages: if both were fed exactly the same signal, zero would result. Can reduce the noise from long leads.

digital. Varying discretely. A digital signal changes to discrete values rather than varying continuously. A digital time display might read 2:57 instead of using the positions of a pair of hands on a clock face.

DIN (Deutsche Industrie Norm). A sort of cable or connector; there are various sorts with different numbers of pins.

envelope form. The overall shape of a signal, outlined by the minimum and maximum recorded values. Often used to display quickly changing signals.

excitation voltage. The voltage supplied to a bridge circuit from which the transducer output signal is derived. Manipulating the transducer changes the measurement elements of the bridge circuit, producing a change in its output voltage.

external trigger. Refers to the connector on the front of the PowerLab marked 'Trigger'. This input lets recording be started from an external source. The trigger level (the voltage necessary to have an effect) depends on the hardware and cannot be changed.

filter. An electronic device or a program that alters data in accordance with specific criteria. Filters in the PowerLab can be used to reduce or to eliminate electronic noise or drift from data readings.

frequency. The number of complete cycles per second of a waveform. Frequency is usually expressed in hertz, Hz (cycles per second), kilohertz, kHz (thousands of cycles per second), or megahertz, MHz (millions of cycles per second).

frequency response. The bandwidth in which a circuit passes a signal without too much attenuation. A low-pass filter's frequency response is the frequency where the output voltage becomes 0.707 ($1/\sqrt{2}$) of the input voltage or has been attenuated by 3 decibels. If a low-pass filter has a frequency response of 200 Hz, say, then the signal is effectively unattenuated up to 150 Hz, and is 0.707 of the original value at 200 Hz.

front-end. An ancillary device that extends PowerLab capabilities, providing additional signal conditioning and features for specialised work. Front-ends are recognised automatically by the PowerLab system and seamlessly integrated into its applications, operating under full software control.

gain. The amount of amplification of a signal.

half-bridge transducer. A bridge transducer only using half of the full-bridge circuit. It consists of two elements of equal value with an excitation voltage applied across them. The output of the transducer is taken at the junction of the two elements.

hertz (Hz). The unit of frequency of vibration or oscillation, defined as the number of cycles per second. The PowerLab's maximum sampling rate can be up to 100 kHz for E series and 200 kHz for SP and ST PowerLabs.

high-pass filter (HPF). A filter that passes high-frequency signals, but filters low ones, by blocking DC voltages and attenuating frequencies below a certain frequency, called the cutoff, or -3 dB, frequency.

I²C. The I²C (eye-squared-sea) connection is used by the PowerLab to control front-ends. It provides power and communications using a 4-wire serial bus (two wires for standard I²C and two control lines).

IEC. International Electrotechnical Commission.

low-pass filter (LPF). A filter that passes low-frequency signals and DC voltages, but filters high ones, by attenuating frequencies above a certain frequency, called the cutoff, or –3 dB, frequency.

Macintosh. A family of Apple computers with built-in graphics and an elegant user interface. Of the Mac OS (operating system) versions, System 7 is needed and System 7.6 recommended normally; a Power Macintosh and Mac OS 8.5 or later are needed for USB.

MacLab. An earlier name for the PowerLab, before it became cross-platform.

Peaks. A simple chromatography analysis package. It reads acquired data from Macintosh Chart, Scope, or text files. (Available for Macintosh only.)

Pod connector. A special 8-pin DIN connector on some PowerLabs giving differential or single-sided connections for some analog inputs (Inputs 1 to 4 on the PowerLab/4SP). Pods can connect to them, and they can also provide power and control for some types of transducers.

Pods. Small, low-cost units that connect to the Pod connectors on the PowerLab. They give alternatives to front-ends for specific tasks, for use with precalibrated transducers and so on.

port. A socket in a computer where you plug in a cable for connection to a network or a peripheral device. Also, any connection for transferring data, for instance between the CPU and main memory.

PowerChrom. A versatile, flexible, and powerful chromatography program for use with a PowerChrom recording unit. (Available for Macintosh only.)

PowerChrom System. The system consists of a PowerChrom recording unit and the

PowerChrom program. It provides generalpurpose manual or automatic recording and analysis of chromatographic peaks.

PowerLab. The PowerLab hardware unit is a self-contained data acquisition hardware unit that connects to a Windows or Macintosh computer. When used in conjunction with programs such as Chart and Scope, it functions as a versatile laboratory instrument.

PowerLab system. The system consists of a hardware unit and applications software (and possibly ancillary devices). It provides a multi-purpose data recording, display, and analysis environment for experimental data.

range. In Chart and Scope, the range is the greatest positive and negative voltage that can be displayed, usually from $\pm 2 \text{ mV}$ to $\pm 10 \text{ V}$, in 12 steps. (Range is inversely proportional to gain, the extent of amplification.)

Scope. An application supplied with a PowerLab that emulates a two-channel storage oscilloscope, with added powerful options. (Macintosh and Windows versions are very similar.)

SCSI (small computer system interface). A connection protocol that provides a computer with fast access to peripheral devices, but has to be set up carefully. SCSI devices can be linked together in a chain, and require termination to prevent signal loss.

serial. A connection protocol for sending information sequentially, one bit at a time, over a single wire. (The PowerLab/800 has a slow serial connection, for use with early Macintosh computers.) **TChart**. A simplified version of the Chart program once used for teaching. Its manual of experiments is now available for Chart.

terminator. A device used in a SCSI chain to maintain signal integrity. Some models of PowerLab have them built in.

transducer. A physical device that converts a mechanical, thermal, or electrical stimulus into a proportional electrical output. For example, there are common transducers to measure force, displacement, temperature, pressure, and similar parameters.


trigger. A signal such as a voltage pulse, used to determine when sampling will begin. Sampling can be made to begin when the trigger level is reached, after it, or even prior to it. See also external trigger.

TTL (transistor-transistor logic). A family of integrated circuits with bipolar circuit logic, used in computers and related devices. TTL is also a standard for interconnecting such ICs, defining the voltages used to represent logical zeroes and ones (binary 0 and 1).

USB (universal serial bus). A relatively new low-to-medium speed connection protocol, for linking multiple peripherals from mice to printers. (Available on new PowerLabs, such as the SP series PowerLabs.)

waveform. The shape of a wave; a graph of a wave's amplitude over time.

Windows. An operating system for PCs with a graphical user interface. Chart for Windows and Scope for Windows require Windows 95 or NT 4 or later. Windows 98, Me, 2000 or later is needed for USB.

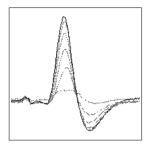

Index

A

ADC (analog-to-digital converter) 28 ADInstruments contacts ii analog input 7–8 analog output 8, 30, 34–35 analog-to-digital converter 29 ASPI 4, 43

B

back panel 9–13 Body protected symbol vii

С

calibration 33 Cardiac protected symbol vii changing fuses 12 Chart 4–5 checking the PowerLab 2 computer requirements Macintosh 4 Windows 3 connections ground 11 I2C 31, 36 Macintosh 23-24 Pod 7-8, 37-38 portable computers 24 PowerBook 23 PowerLab to computer 17-26 safe, to humans 5, 7, 31 safety v-viii SCSI 20-26, 36 USB 17-20, 35 Windows 24-26

contacting ADInstruments ii cooling fan 13

D

DAC (digital-to-analog converter) 30 DC drift 32 digital input 31, 37 digital output 31, 37 digital-to-analog converter 30

E

equipotential connection 11 error patterns 16–17 external trigger 9, 29–30, 33

F

fan 13 FIFO devices 27, 29 front panel 5–9 front-ends 5 fuse holder 12 fuses, changing 12

G

ground connection 11

I

I2C port 8, 10, 36, 38 installation problems 43

M

Macintosh PowerBook 23 maintenance viii

N

New Hardware wizard 43

P

Pod connection 7-8, 37-38 Pods 8 portable computers 24 Power indicator 7 power switch 12 PowerLab accuracy 33 back panel 9-13 calibration 33 connectors 35-38 front panel 5–9 input amplifiers 31 power supply 31 safety v-viii, 5, 7, 31 self-test 16-17 user modification 27 voltage rating 3 PowerLab system 3 problems and solutions Macintosh 39-42 Windows 42–46

R

replacing fuses 12

S

safety v—viii, 5, 7, 31 Scope 4–5 SCSI cables 21, 22, 23 card 23–26, 43 chain 22 connection 20–26, 36 connection rules ??–20, 20–22 ID number 21, 22 multiple devices 24, 25 port 10, 35 problems 39–46 termination rules 23 SCSI ID switch 10, 22 SCSIProbe 22 self-test 16–17 Status indicator 7, 23 storage vii

T

technical specifications PowerLab/16sp 53–56 PowerLab/4sp 47–50 PowerLab/8sp 50–53 termination rules for SCSI 23 transistor-transistor logic 33 trigger 9 Trigger indicator 9 TTL (transistor-transistor logic) 33

U


USB cables 20 connection 17–20, 35 connection rules 19–20 port 9 user modification 27

V

vent holes 13 voltage rating 2–3

W

Warning Symbol vii

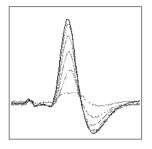

Licensing & Warranty Agreement

Extent

This Agreement is between ADInstruments Pty Ltd ['ADI'] and the purchaser ['the Purchaser'] of any ADI product — software, hardware, or both and covers all obligations and liabilities on the part of ADI, the Purchaser, and other users of the product. The Purchaser (or any user) accepts the terms of this Agreement by using the product. Any changes to this Agreement must be recorded in writing and have ADI's and the Purchaser's consent.

Copyright and Trademarks

ADI develops proprietary computer software and hardware, including MacLab[®], PowerChrom[®], and PowerLab[®] units. All ADI software, hardware, and documentation is protected by copyright, and may not be reproduced or altered in any way, nor may products be derived from or based on it. ADI retains the exclusive ownership of the trademarks and registered trademarks represented by its company name, logo, and product names.

Responsibilities

The Purchaser and any others using any ADI product agree to use it in a sensible manner for purposes for which it is suited, and agree to take responsibility for their actions and the results of their actions. If problems arise with an ADI product, ADI will make all reasonable efforts to fix them. This service may incur a charge, depending on the nature of the problems, and is subject to the other conditions in this Agreement.

Limitations

ADI products are produced to high standards, and should perform substantially as described in the supplied documentation. There is a limited hardware warranty, and technical support is provided for all products.

Nevertheless, since ADI products could be affected by external factors (for instance, the computer system on which they run), absolute performance and reliability cannot be guaranteed. No warranty, either express or implied or statutory, other than that contained in this Agreement, is made in respect to ADI products. The Purchaser therefore assumes all risks as to the performance and reliability of the products, and the results gained using them. ADI is not responsible for any problems with the computer system not directly related to ADI products.

No claim for damages of any sort arising from the use or misuse of ADI products can be made against ADI or its agents or employees.

Hardware Warranty

ADI warrants that the hardware purchased by the Purchaser shall be free of defects in material and workmanship for one year from its date of purchase. If there is such a defect, ADI will repair or replace the equipment as appropriate, and the duration of the warranty shall be extended by the length of time needed for repair or replacement.

To obtain service under this warranty, the Purchaser must notify ADI of the defect before the warranty expires. Unless other arrangements are made, ADI will provide an RMA (Return Merchandise Authority) number, and a service centre address to which the Purchaser must ship the defective product at his or her own expense. The product should be packed safely (preferably in its original packaging) and have the RMA number on the shipping label, beneath the address. ADI will pay return shipping.

This warranty applies only to hardware used normally and within its operating specifications. It does not cover hardware modified in any way, subjected to unusual physical, electrical, or environmental stress, used with incorrectly wired or substandard connectors or cables, or with the original identification marks altered.

Software License

The Purchaser has the non-exclusive right to use the supplied ADI software. (The Purchaser's employees or students, for instance, are entitled to use it, provided they adhere to this Agreement.)

The Purchaser can make as many backup copies of ADI software as is reasonable. Each separate purchase of a software program, however, licenses it to be used on only one computer at any given time. Although multiple copies of a purchased program may exist, two or more copies must not be used simultaneously. Site licenses (multipleuser licenses) are treated as if five copies of a program were purchased separately, even if only one CD or set of disks is provided.

Technical Support

Once a customer registration form has been filled out and sent in, the Purchaser is entitled to free technical support for any ADI product for one year from its date of purchase. (A customer registration form should be included with each product: if for some reason you cannot find one, contact ADI for another.) This involves help and advice concerning installation, operation, particular uses, and problems with ADI products. Services not directly related to ADI products may incur a charge.

Jurisdiction

This Agreement shall be governed by the laws of New South Wales in Australia, and any proceedings concerning it shall be heard and resolved by the Supreme Court of New South Wales in Australia.